



**replacement parts industries, inc.**

*"The Alternate Source"®*

CALL 800-221-9723 or 818-882-8611 • FAX 818-882-7028

E-MAIL: [techsupport@rpiparts.com](mailto:techsupport@rpiparts.com) • WEBSITE: [www.rpiparts.com](http://www.rpiparts.com)

## **CAC016 SPEED CONTROL ADJUSTMENT INSTRUCTIONS**

1. Install speed control in normal manner.
2. Turn shaft on potentiometer counter-clockwise until it stops.
3. Press the enclosed knob onto the potentiometer shaft.
4. Included with the knob is an "indicator line" with adhesive on one side. Locate the "zero" position on the front panel of the centrifuge. Remove the "indicator line" from back-up paper and press onto the knob, making sure it points to zero.