



**replacement parts industries, inc.**

*"The Alternate Source"®*

CALL 800-221-9723 or 818-882-8611 • FAX 818-882-7028

E-MAIL: techsupport@rpiparts.com • WEBSITE: www.rpiparts.com

## **IEC016**

*(Old RPI Part #DAC016)*

# **RUBBER CUSHION INSTALLATION INSTRUCTIONS**

*(Must be used if micro-capillary hematocrit tubes are sealed in plasticene.)*

1. The inner side of the flange of the head should be lined with the cushion (RPI Part #IEC016). The ends should make a smooth butt joint (cut as necessary to obtain correct length). The butt joint should not be opposite a groove in the head.
2. As the capillary tubes begin puncturing the cushion, move it so that there is always ample cushion material behind the tubes. Replace entire cushion when this is no longer possible.